



APPLICATIONS

- CIP/SIP System Condensate Drainage
- Sterilization of Process Vessels
- Culinary Steam
- Humidifiers
- WFI System Sterilization
- Fermenter Sterilization

OPTIONS

- MP Mechanical Polish to 10 μ in. (0.25 μm) Ra
- EP Electropolish
- SLR SLR Orifice
- Tef-Steel, PTFE, Teflon®, E.P.D.M., USP Class VI & other gasket materials available
- -B Bellows for low subcool
- 6" Condensate Leg (CME-W25, CME-A)

APPLICABLE CODES

ASME BPE

Canadian Registration # 0E0591.9

CDS SANITARY **THERMOSTATIC STEAM TRAPS**

Pressures to 100 PSIG (6.9 barg) Temperatures to 338°F (170°C)

Steepest Interior Surfaces—Designed to completely drain without puddling.

Stainless Steel Body—Body Material is 316L Stainless Steel with 20 μ in. (0.5 μ m) Ra internal finish and 32 μ in. (0.75 µm) Ra external finish. Available with mechanical polishing to 10 µ in. (0.25 µm) Ra and/or electropolish.

Self centering Valve—Leak tight shut off. Assembly of actuator and valve to impingement plate allows the valve to self align with center of the orifice.

Temperature Sensitive Actuator—One moving part. 316L Stainless Steel, fail open, welded actuator for maximum corrosion, thermal and hydraulic shock resistance.

Directional Discharge—Erosion prevented by directing discharge to center of piping.

Maintenance—Can be easily removed and disassembled for sterilization and/or repair.

Three Year Guarantee—Guaranteed for three years against defects in material or workmanship.

Industry Standard Food Grade Gasket—White Viton food grade gasket offers superior performance for higher pressure steam applications.

Large Orifice Selection—Broad selection of orifice sizes provide greatest sizing and selection flexibility.

Superior Air Handling—Best air handling capability provides for fast startup.

Unique SLR Orifice OptionSpecify when immediate elimination of condensate and improved sensitivity is desired. An orifice on the valve allows for continuous discharge of condensate. Trap will nominally pass 50 lb/hr (22.7 kg/hr) of condensate at 50 psi (3.4 barg) within 0.5°C (1°F) of saturated temperature.

Bar Stock—Connectiion fittings are not welded onto inlet and outlet pieces.

Models

- CDS202-Low capacity
- CDS203–Medium capacity
- CDS204-High capacity

NOTE: Please specify if Material Test Reports (MTR) or Certificates of Conformance (COC) are required.

OPERATION

Thermal actuator is filled at its free length with a liquid having a lower boiling point than water. On start-up, valve is normally open to discharge air, non-condensibles and condensate. When steam enters trap, thermal actuator fill vaporizes to a pressure higher than line pressure. This forces

valve into seat orifice to prevent any further flow. As condensate collects, it takes heat from the actuator, lowering internal pressure. Line pressure will then compress thermal actuator to open valve and discharge condensate. Valve opening automatically adjusts to load conditions from minimum on very light loads to full lift at maximum load.

CDS SANITARY THERMOSTATIC **STEAM TRAPS**

SPECIFICATION

Steam trap shall be of balanced pressure design with 316L welded bellows capable of releasing condensate within 1°C (2-3°F) or 5°C (10°F) of saturated pressure. All other interior wetted components shall be of 316L stainless. It shall have interior body finish of at least 20 μ in. (0.5 μm) Ra and exterior body finish of at least 32 μ in. (0.75 μm) Ra. Trap shall utilize hygienic body clamp allowing disassembly for inspection or cleaning and be entirely self draining when installed in vertical configuration. Trap end connections shall be standard hygienic clamp. Thermostatic actuator shall employ a conical valve lapped to the seat. A minimum of three orifices shall be available. Traps shall have SLR orifice where drainage at saturated temperatures is required. Traps shall be guaranteed against defects for 3 years.

MAXIMUM OPERATING CONDITIONS

PMO: Max. Operating Pressure 100 psig (6.9 barg) TMO: Max. Operating Temperature 338°F (170°C) 150 psig (10.3 barg) PMA: Max. Allowable Pressure TMA: Max. Allowable Temperature 366°F (186°C)

BODY SURFACE FINISH

Internal <20 μ in. (0.5 μ m) Ra SFCI.External <32 μ in. (0.75 μ m) Ra. Optional mechanical polishing to 10 µ in. (0.25 µm) Ra and/or electropolish SFC4

GASKET APPROVALS

FDA CFR Title 21 Part 177, USDA, USP Class VI, 3A Sanitary Standard, NSF

SERVICE NOTES

Trap is designed to be self draining for vertical installation (discharge down) 1/2" - 3/4" service trap should be installed with 3/4" inlet gasket. 1" - 11/2" service trap should be installed with 11/2" inlet gasket.

SLR ORIFICE OPTION


Specify when immediate elimination of condensate and improved sensitivity is desired. An orifice on the valve allows for continuous discharge of condensate. Trap will nominally pass 50 lb/hr (22.7 kg/hr) of condensate at 50 psi (3.4 barg) within 0.5°C (1°F) of saturated temperature.

Polishing Procedure

All surface finishes are achieved without the use of additional buffing, compounds or grit.

-B Bellows

1°C (2-3°F) subcool for sensitive applications under 40 psi (2.7 barg)

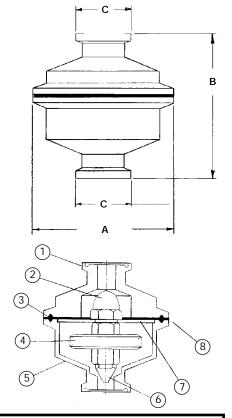

Connections: 1/2" - 11/2" Hygienic Clamp

DIMENSIONS inches (mm) and WEIGHTS pounds (kg)

Service	Α	В	С	Weight
1/ on al 3/	2½	2%	63/64	1.8
½ and ¾	(63)	(67)	(25)	(8.)
1 and 11/	2½	25/8	163/64	2.3
1 and 1½	(63)	(67)	(50)	(1.0)

MATERIALS OF CONSTRUCTION				
Item	Part Name	Material		
1	Body - Inlet	316L A276		
2	Actuator Nut	316L		
3	Gasket	Viton 3227		
4	Actuator	316L		
5	Body - Outlet	316L A276		
6	Valve	316L		
7	Clamp (not shown)	304		
8	Impingement Plate	316L		

Maximum Capacity—lbs/hr 10°F Below Saturation (Kg/hr 5°C Below Saturation) Differential PSIG (bar) Orifice Trap 30 40 50 60 70 80 100 10 20 (0.34)Inches (0.7)(1.4)(2.1)(2.8)(3.4)(4.2)(4.9)(5.6)(6.2)(6.9)291 411 581 719 831 919 1000 1075 1130 1174 1207 CDS 202 5/32 (377)(132)(186)(454)(488)(533)(547)(264)(326)(417)(513)825 1495 1975 2175 2825 550 1210 1750 2350 2525 2650 CDS 203 1/4 (249)(374)(549)(678)(794)(896)(987)(1066)(1145)(1202)(1281)1671 2128 2301 2509 2713 2819 2946 3091 3159 1520 1869 CDS 204 5/16 (689)(758)(848)(965)(1044)(1138)(1230)(1278)(1336)(1402)(1433)

NOTE: Please specify if Material Test Reports (MTR) or Certificates of Conformance (COC) are required.